


Curie Point Injector JCI-22

Odor Analysis Fried rice cracker



Air Sampler
JAS-15M II

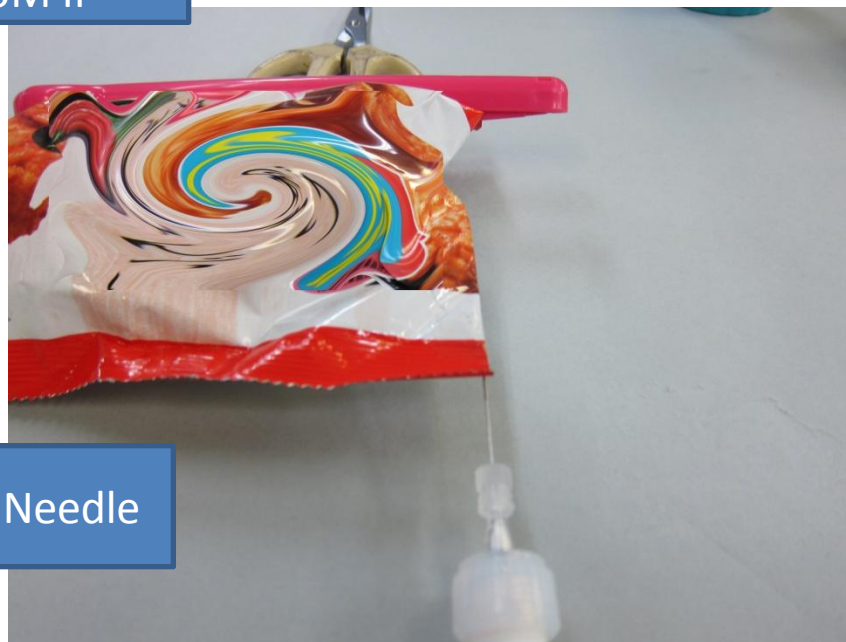

Sampling Needle

Trap tube
mini-PAT



JCI-22
JAS-15M II

Japan Analytical Industry Co., Ltd


Odor Analysis

Fried rice cracker



mini-PAT

Quartz Glass Sample Tube
ID: 4mm Length: 36mm
Absorbent :Tenax GR 25mg
Pyrofoil :F280 (280c)

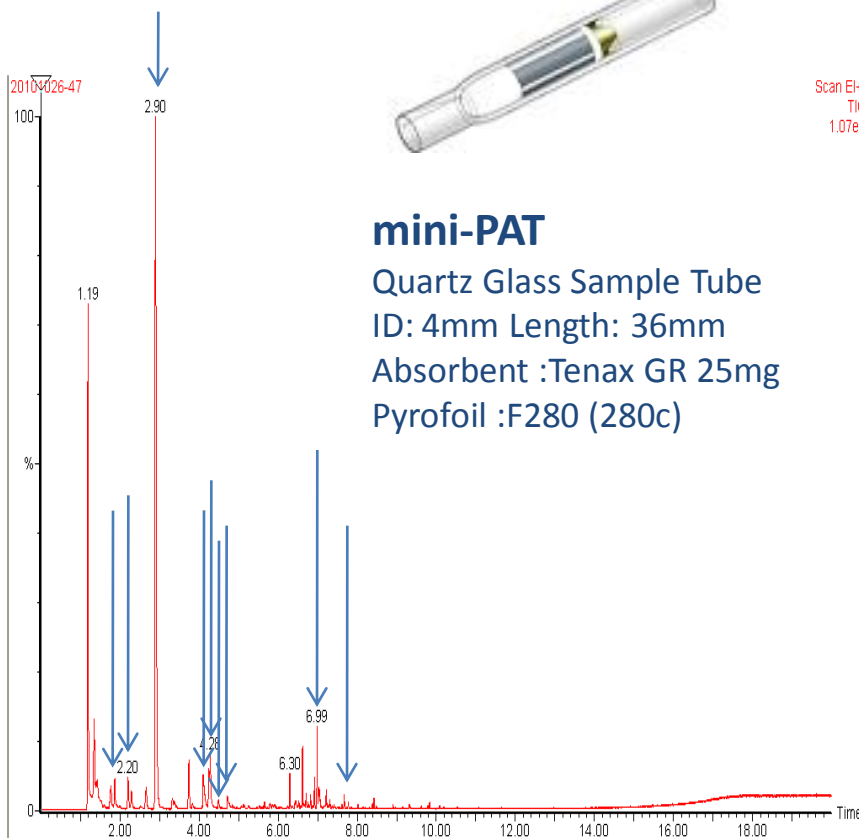

Clarus 600T GC/MS

- Anal time : 20min
- Oven temp : 40°C(3min)-(20°C/min)-320°C(3min)

MS Condition

- Ion source Temp : 250°C
- Interface Temp : 300°C

Column

- DB-5 MS: 30 m, 0.25 mm I.D, 0.25 um df.
- TIC

From early retention time →

Methylbutylketone

Isovaleraldehyde

Propylacetate

Butanediol

Hexanal

Ethylacetate

Purimidinemethyl

Ethylhexanol

Phenylethanol



JCI-22
JAS-15M II

Japan Analytical Industry Co., Ltd